Compact Dry ETB medium for enterobacteriaceae/Enterobacteriaceae/milieu pour nombre enterobacteriaca selectifs /medio para enterobacterias/ medio per conta enterobacterias /meio para contagem enterobacterias

40 plates/Platten/plaques/placas/piastre/placas

ID-No. 1 002 941

240 plates /Platten/plaques/placas/piastre/placas

ID-No. 1 002 942

920 plates/Platten/plaques/placas/piastre/placas

ID-No. 1 002 943

English

Deutsch

Français

Compact Dry ETB is a ready to use, selective plate containing glucose for the enumeration of Enterobacteriaceae

Specimen pretreatment

Viable count in water or liquid foodstuff

Drop 1 ml of specimen (dilute if necessary) on the middle of the Compact Dry plate.

Viable count in solid foodstuff

Add buffer solution to the specimen and homogenize by stomacher®. Drop 1 ml of specimen (dilute if necessary) on the middle of the dry sheet of the Compact Dry plate.

Viable count in swab test specimen

Use the swab to wipe the surface, put into the device with wiping solution. Drop 1 ml of wiping solution (dilute if necessary) on the middle of the Compact Dry plate. It is recommended to use "Swab for Compact Dry" offered by HyServe Id-No. 1 002 952/3 (40/240 pieces).

Test instructions

1. Open the cap and drop 1 ml of specimen on the middle of the Compact Dry plate.
2. Specimen diffuses automatically and evenly into the sheet and transforms the dried sheet into a gel within seconds.
3. Put the cap again on the plate and write the information needed on the memorandum section.
4. Turn over the capped plate and put in the incubator.
5. After incubation count the number of **red/purple** colored colonies underneath the plate. White paper placed under the plate helps to count the colonies.

Incubation time 24 ± 2 hours

Incubation temperature 37 ± 1°C

Interpretation of the results

Colonies grown are almost all **red/purple**. **Red/purple** colored colonies are Enterobacteriaceae. **Red/purple** colored colonies and otherwise colored colonies together are the total gram negative bacteria.

Storage and shelf life

Keep at room temperature (+ 1 to +30 °C).

Total shelf life 16 months after manufacturing.

Notes

- Some colonies might not be clearly **Red/purple** colored.
- High concentrations on plates will cause the entire growth area to become red/pink. In this case dilute the specimen.
- After use please follow the current disposal regulations.
- The growth area is 20 cm². The back of the plate has a grid carved of 1 cm x 1 cm to make the colony counting easier. In case of any difficulties to count colonies due to large number of colonies grown, the bacteria count can be obtained by multiplying 20 by an average number of colonies per grid counted from several grids. However, dilutions are recommended.
- Compact Dry plates are produced at an ISO 9001 /ISO 13485: 2003 certified site.
- MicroVal approval No.0806-006 LR according to ISO 21528 (2004) and ISO EN 16140:2003
- NordVal approval No. 034

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